Reinigungshinweise miniPERM®

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Cleaning instructions miniPERM®

Please read the following information before using the miniPERM®!

⚠ For disinfection, use only 70 % pharmaceutical grade ethanol or Isopropanol for the miniPERM® bioreactor, the Universal Turning Device, the incubator and water bath.

Do not use n-propanol!

⚠ All other chemicals, especially quarternary ammonium complexes, may cause microcracks.

Cleaning instructions for the autoclavable nutrient module and the autoclavable screw cap of the miniPERM®

Autoclavable nutrient module

- Clean only with hot water, and neutral or weak acid or alkaline detergent (e.g. RBS Neutral, Roth GmbH) without rinsing agents.
- Rinse with ample water to completely remove the detergent.

⚠ Do not use an automatic dish washer!

- Autoclave without screw cap to 121°C for 20 min (max. 10 times).

⚠ Do not use any alkaline-based corrosion protector in autoclave boiler water!

Autoclavable screw cap with integrated gas exchange membrane

- Rinse screw cap with distilled water only.
- Air dry and wrap in appropriate bag or film prior to autoclaving.
- Autoclave to 121°C for 20 min (max. 5 times).

⚠ Do not use ethanol, propanol or disinfectants / solvents that contain alcohol, as their use will impair the function of the PTFE membrane!